



**BLDEA's SSM College of Pharmacy and Research Centre**  
**Vijayapur - 586103**  
**Dept. of Pharmacology**

**STANDARD OPERATING PROCEDURE**  
Updated: 01-02-2022 Valid upto 01-08-2022

# HOTAIROVEN

## Standard Operating Procedure

- Load the properly packed material to be sterilized within the oven,
- Close the door.
- Connect the plug to the power source and switch on the power.
- Switch on the mains of oven.
- LED lamp for the power indicator will glow
- Set required temperature (160oC for 2 hours for sterilization of glassware) using temperature adjustment knobs.
- Heaters will start, the LED for heating will start glowing and the temperature rises steadily to the Set temperature.
- Note the time at which the set temperature is reached.
- After completion of sterilization, switch off the MAINS and the power source.

**Prepared by: Shri Shridharkumar.S. Biradar**

**Checked by: Dr. V P Patil**

**Approved by: Dr. Shivakumar Hugar**