


**STANDARD OPERATING PROCEDURE OF**  
**MOTAR STIRRER**

 <p><b>BLDEA'S Shri Sanganabasava Mahaswamiji College of Pharmacy &amp; Research Centre Vijayapur 586103</b></p>	<b>DEPARTMENT: PHARMACEUTICS</b>	<b>SOP NO:</b>
	<b>AREA: PHARMACEUTICS LABORATORY</b>	<b>Company Name: REMI lab Model no: INSUB</b>
	<b>SUBJECT: SOP FOR MOTAR STIRRER APPARATUS</b>	<b>PAGE NO: 1-2</b>
		<b>EFFECTIVE DATE: 25/02/2022</b>
		<b>REVIEW PERIOD: 25/02/2023</b>

**Purpose:**

To provide a procedure for the operating of Motor stirrer apparatus

**Scope:**

Applicable to operation of Motor stirrer apparatus

**References:**

Equipment Manual

**Safety Issues and Precautions:**

- Wore Head cap, Mask, Hand gloves and other safety requirements during the performance of stated activity
- Please make the entry of Usage with require details in the Equipment log book
- Handle the Instrument properly with care

**Procedure:**

- 1) Production person shall ensure the cleanliness of the stirrer
- 2) Fix the shaft or propeller to the mortar. Keep the shaft into the suspension to the half of the depth of suspension or up to that level so that suspension should not flush out
- 3) Put ON the electrical supply and start the stirring

- 4) Stirring is carried out until a uniform slurry is obtained or time specified in BMR
- 5) Switch OFF the mains after completion of the operation
- 6) At fix to be cleaned label and fill the equipment log book

	<b>PREPARED BY</b>	<b>CHECKED BY</b>	<b>APPROVED BY</b>
<b>NAME</b>	Mrs. Ashwini S. Gudigenavar	Mr. A B. Walikar	Dr. S C. Marapur
<b>DESIGNATION</b>	Asst. Professor	Asst. Professor	Professor & Head
<b>DEPARTMENT</b>	Pharmaceutics	Pharmaceutics	Pharmaceutics
<b>SIGNATURE &amp; DATE</b>			

