



BLDEA's SSM College of Pharmacy and Research Centre
Vijayapur - 586103
Dept. of Pharmacology

STANDARD OPERATING PROCEDURE
Updated: 01-02-2022 Valid upto 01-08-2022

POLE CLIMBING APPARATUS

FUNCTION: To study the anti psychotic activity by using pole climbing apparatus

Standard Operating Procedure:

- Connect the power connection and switch on the instrument.
- Adjust the knob for sound or light.
- Time for buzzers or light.
- Adjust voltage of shock to be given.
- Open upper lid and animal on the grid floor for the observations.
- Push the buzzers or light on a sound or light will come for specified time and response.
- Now apply shock and note response.
- Repeat this for several times.
- Switch Off the main and remove animals.

Prepared by: Dr.H.M.Nanjappiah

Checked by: Mr. V P Patil

Approved by: Dr. Shivakumar Hugar