



BLDEA's SSM College of Pharmacy and Research Centre
Vijayapur - 586103
Dept. of Pharmacology

STANDARD OPERATING PROCEDURE
Updated: 01-02-2022 Valid upto 01-08-2022

DIGITAL TELE THERMOMETER

FUNCTION: To study the antipyretic effect.

- **Standard Operating Procedure**
- Connect the rectal probes into their respective sockets of the telethermometer by matching their numbers.
- Insert the probe into the rectum of the animals to record rectal temperature of the control and test group challenged with Baker's yeast.
- Allow sufficient time to attain equilibrium so that the reading gets stabilized.
- Note the rectal temperature of the animal individually.

Prepared by: Shri Shridharkumar S. Biradar

Checked by: Dr. V.P.Patil

Approved by: Dr. Shivakumar Hugar