



**BLDEA's SSM College of Pharmacy and Research Centre**  
**Vijayapur - 586103**  
**Dept. of Pharmacology**

## **STANDARD OPERATING PROCEDURE**

Updated: 01-02-2022 Valid upto 01-08-2022

# **RADIAL ARM MAZE**

### **Standard Operating Procedure**

1. Train rats in radial maze.
2. Keep food pellets at the end of arm
3. Place the rat at the centre of the maze
4. Record observations (Time taken by rat to find food pellet, frequency of defecation and urination)
5. Allow the rat to explore maze for 5-10 minutes as per study taken
6. Perform trial for all rats of Control and test group

**Prepared by: Shri Shridharkumar. S. Biradar**

**Checked by: Dr. V P Patil**

**Approved by: Dr. Shivakumar Hugar**